NICHIREI

N-Histofine® **Simple Stain AEC Solution** Peroxidase Chromogen/Substrate Solution

500 tests **Code: 415182F**
1500 tests **Code: 415184F**

AMINOETHYL CARBAZOLE (AEC) SUBSTRATE SOLUTION **SUBSTRATE FOR HORSERADISH PEROXIDASE**

I-1. DESCRIPTION

N-Histofine® Simple Stain AEC Solution is designed as one of reagents to visualize antigen binding primary antibody on tissue sections.

I-2. INTENDED USE

N-Histofine® Simple Stain AEC Solution is designed to allow immunohistochemical tests, to reveal antigens that reacts with a user-supplied primary on tissues and cells. The product is characterized by a coloring reagent/substrate for coloring a detection system reagent labeled with peroxidase. The product is characterized by a coloring reagent/substrate for coloring a detection system reagent labeled with peroxidase. The product is one of the specific chromogen/substrate solutions for peroxidase and develops red colored precipitate by reaction with peroxidase. The result provides a physiological or pathological state.

For In-vitro Diagnostic Use. For manual staining. Staining Results are qualitative. For formalin-fixed paraffin-embedded human tissue sections.

For professional users.

Interpretation must be made within the context of the patient's clinical history and other diagnostic tests by a qualified pathologist.

I-3. PRINCIPLE

The antigen / antibody / Universal Immuno-peroxidase Polymer complex can be prepared by allowing the reagent to react with a primary antibody bound to the antigen on tissue section. The enzymatic activity of this complex results in a colored deposit, thus staining the antigen site.

II. COMPONENTS

Liquid. Ready to use.

Simple Stain AEC Solution	[500 tests]	[1500 tests]
...	17 ml × 3 bottles	17 ml × 9 bottles

III. USAGE

Adjust **N-Histofine®** Simple Stain AEC Solution to room temperature prior to use.

React the objective tissue section with peroxidase conjugate in advance.

At the same time, staining of Negative control slide using PBS in place of primary antibody is recommended to be prepared for checking the staining result.

For color development, incubate the tissue section on the slide with 1-2 drops of **N-Histofine®** Simple Stain AEC Solution using Dropping bottle at room temperature for 5-20 minutes.

Adjust the reaction time by microscopic observation by contrast with the Negative control slide.

After adequate coloration, wash the slide with tap water.

This red deposit is soluble in organic solvent. Therefore, after counterstain, the slide can be mounted with water-soluble mounting media.

IV. STAINING PROCEDURE

Refer to the instruction* of **N-Histofine®** Simple Stain MAX PO (MULTI) for staining procedures such as Principle, Materials required but not supplied, Specimen preparation, Control slides, Interpretation of results and General limitation. Validation is required when other detection systems besides above are used.

*Download of the instruction is available at following website:

http://www.nichirei.co.jp/bio/english/products/pdf/CE_MAXPO_MULTI.pdf

V. STORAGE, HANDLING & SHELF LIFE

1. Store at 2-8°C.
2. Confirm expiration date indicated on the labels of vial and carton before use.
3. Adjust this product to room temperature before use.
4. Refrigerate this product promptly after use.
5. The reagent is stable 18 months after manufacturing.

VI. PRECAUTIONS

1. Before using this reagent, read these instructions.
2. Do not use reagents after the expiration date.
3. For professional users.
4. Specimens, before and after fixation, and all other materials exposed to them, should be handled like biohazardous materials with proper precautions.
5. Inhalation or ingestion of the highly allergic fixative formaldehyde is harmful. Wear protective mask. If swallowed, induce vomiting. If skin or eye contact occurs, wash thoroughly with water.
6. Organic reagents are flammable. Do not use near open flame.
7. Never pipette reagents by mouth and avoid their contact with skin, mucous membranes and clothes.
8. Avoid microbial contamination of reagents as incorrect result may occur.
9. Avoid splashing of reagents or generation of aerosols.
10. Unused solution should be disposed of according to local, State and Federal regulations.
11. Any serious incident that has occurred in relation to **N**-Histofine® Simple Stain AEC Solution shall be reported to the manufacturer and the competent authority of the Member State in which the user and/or the patient is established.

VII. REFERENCE

- (1) Vilayvong S, et al. Useful Application of Immunostaining to Malignant Pleural Effusion among Lao People in Vientiane Capital, Lao PDR. Asian Pac J Cancer Prev. 2019 Jan 25;20(1):243-248.
- (2) Rochette A, et al. Asporin is a stromally expressed marker associated with prostate cancer progression. Br J Cancer. 2017 Mar 14;116(6):775-784.
- (3) Saho S, et al. Active Secretion of Dimerized S100A11 Induced by the Peroxisome in Mesothelioma Cells. Cancer Microenviron. 2016 Dec;9(2-3):93-105.

 REF	Catalog/Code Number		Temperature Limitations	 IVD	In Vitro Diagnostic Medical Device
	Manufacturer	 LOT	Batch Code		Contains Sufficient for <N> Tests
	Use By		Consult Instructions for Use	 EC REP	Authorized Representative in the European Community
	CE-mark, code of the notified body		For IVD Performance Evaluation only	 SAMPLE	Sample

NICHIREI BIOSCIENCES INC.
6-19-20, Tsukiji, Chuo-ku, Tokyo 104-8402, JAPAN
Phone:81-3-3248-2208, Facsimile:81-3-3248-2243

MedEnvoy Global B.V.
Prinses Margrietplantsoen 33-Suite 123
2595 AM The Hague
The Netherlands

N-Histofine[®]
Simple Stain AEC Solution
 Peroxidase Chromogen/Substrate Solution

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